


Dip Sampler

for liquids

Disposable Plunging Siphon

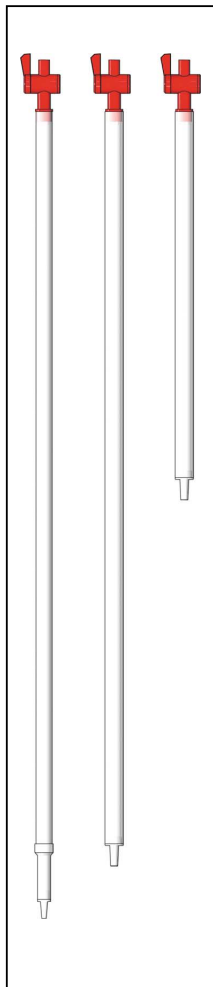

Dip Sampler


made of PP for rapid and simple spot checks of liquids. The dip samplers are transparent and enable the contents to be seen clearly. Upper end with reflux valve, lower end with ball valve.

- Transparent PP
- Type 200 variable step by step from 100 to 200 cm
- Special sizes on request
- All-layer sample and target sample possible

Operation

All-layer sample 1. Insert the dip sampler with **opened valve** slowly into the liquid. 2. Close the valve. 3. Withdraw the dip sampler. Remove the sample.

Target sample 1. Insert the dip sampler with **closed valve** into the liquid until the target spot. 2. Open the valve. The liquid flows in. 3. Close the valve. 4. Withdraw the dip sampler. Remove the sample.

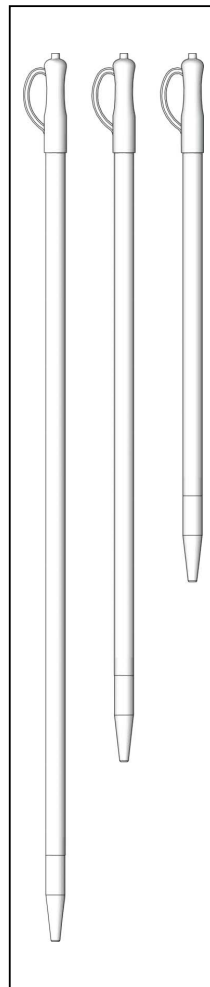

Disposable Plunging Siphon

If strongly adhesive liquids are sampled, it can be a lot less expensive to dispose of the sampler properly than to employ a long and complicate cleaning process.

The disposable plunging siphons were specially developed for this. They are made of PE, transparent and allow easy recognition of their contents. PE provides a surface that resists adhesion of the sample material, minimising sample retention and loss.

Operation

1. Insert/fill collector into liquid.
2. Close thumbhole above the handle using your thumb.
3. Remove plunging siphon, and empty liquid into the sampling bottle.

Dip Sampler



Length cm	Diameter mm	Price/Unit	Price/Unit of (X) units	Article No.
50	22	30,80	(10) 28,10	5615-50 N4
100	22	33,20	(10) 30,20	5615-100 N4
200 ¹⁾	22	39,80	(10) 35,80	5615-200 N4

¹⁾ variable step by step from 100 cm to 200 cm

Disposable Plunging Siphon, complete with handle



Total length cm	Plunging depth cm	Capac- ity ml	Tube diameter mm	Price/ unit	Price/unit of (X) units	Article No.
50	35	75	27	6,20	(25) 5,60	5387-50 N4
75	60	175	27	11,30	(25) 10,10	5387-75 N4
100	85	275	27	16,30	(25) 14,70	5387-100 N4

Prices in € – Phone + 49 (0) 7621 9331-16, Fax + 49 (0) 7621 9331-31