


Zone Sampler

for all kinds of bulk goods, according to DIN 53 242
for fast cross-section samples

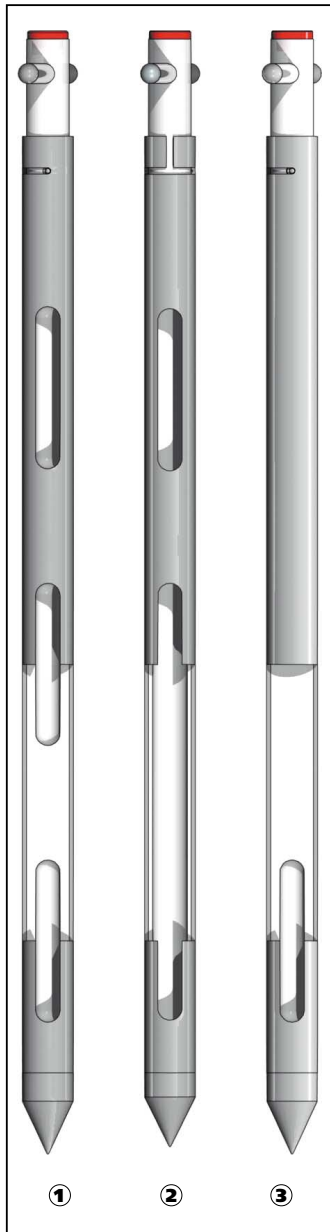

Zone Sampler

Three versions of the Zone sampler are available for different sampling tasks.

1 Multi-sampler with several closed chambers for multi-sampling from several predefined sample depths.

2 All-layer sampler and Jumbo sampler with open inner tube, no chambers, for cross-section samples through all layers of the sample. Simple emptying through open end of sample tube. The Jumbo sampler with a tube diameter of 50 mm is ideal for larger grain sizes and amounts.

3 Uno-sampler with one closed chamber, for a targeted sample from a specific sample depth.



Color coding

The colored push-in buttons (sterilizable) supplied as standard can be used to identify Multi-samplers and Uno-samplers to avoid mix-ups.

Bürkle Zone Samplers

are used for sampling all types of bulk goods. The great advantage of the zone sampler is that **cross-sectional samples** can be taken from all the layers of a container.

The Zone samplers can take samples from a **wide range of bulk goods**, from the finest powders up to coarse grains such as corn or nuts.

The Zone samplers are available in **4 different lengths**. 55 cm - specially suitable for sacks, 85 cm - specially for barrels, drums and mixing containers, and 150 cm and 250 cm for Big-Bags, silos and tanks. Special lengths are available on request.

Safe and reliable sampling is made even easier with useful special accessories such as sealing sleeves, Quali-Paper, cleaning brushes and close-it control seals.

Advantages

- Smallest possible gap between inner and outer tube
- Easy cleaning, removable tip
- Color coding (Multi- and Uno-samplers)
- Ultrapure materials: stainless steel 1.4306/PTFE
- 4 lengths from 55 to 250 cm
- Outstanding finishing, robust execution

Operation

1. Turn handle to close chamber.
2. Push in sampler to required depth.
3. Turn handle to open chamber, sample flows into chamber.
4. Turn handle to close chamber again.
5. Pull out sampler.
6. Empty chamber onto Quali-Paper.
7. Close container opening with close-it (see pages 92ff).

seal-it® Security seal for sample containers

- Security provided by unique 7-digit control number
- Fits any bottle, can, bag, container
- Additional documentation label with same number
- Detailed information on page 105

Test instruments

Would you like to carry out a test before making a purchase? We will be pleased to send you a test instrument on request. A small charge will be made for maintenance and cleaning work.